


# The Perfect Qubit: How Tunable Light Helps in the Quantum Quest

Widely tunable continuous-wave lasers based on OPO technology make it quicker and easier to characterize different qubit contenders

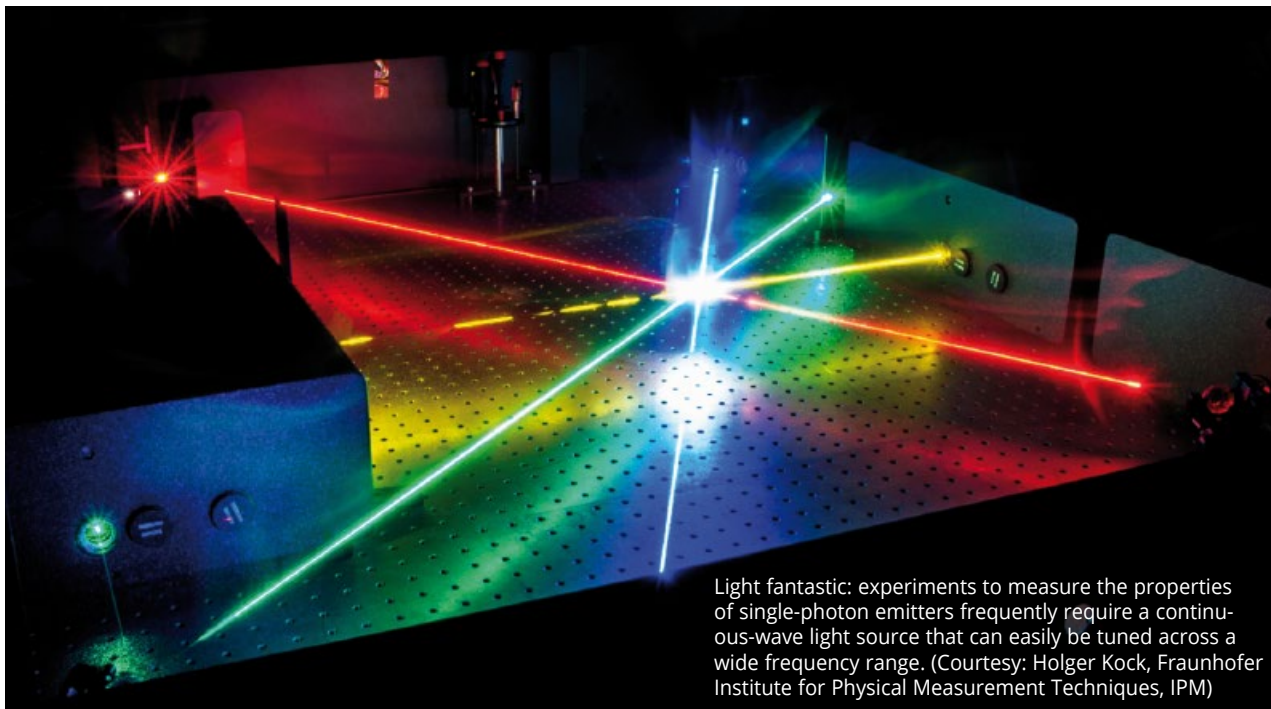

Light fantastic: experiments to measure the properties of single-photon emitters frequently require a continuous-wave light source that can easily be tuned across a wide frequency range. (Courtesy: Holger Kock, Fraunhofer Institute for Physical Measurement Techniques, IPM)

Physicists are still looking for the perfect quantum qubit: a two-level quantum system that can be precisely measured and controlled, while remaining unaffected by its environment. One of the most promising candidates are defect centers in solid-state materials, also known as colour centers, which have been found to emit a single photon per excitation event when excited by laser light of a particular frequency.

Most early attention has focused on nitrogen vacancy centers in diamond – which offer single-photon emission at room temperature – but they are not ideal for all applications because their asymmetric charge distribution makes them sensitive to local fluctuations in the electric field.

Researchers are therefore investigating the properties of different colour centers in diamond, including silicon and germanium vacancies, as well as other material systems such as 2D hexagonal boron nitride (hBN). But it can be difficult and time consuming to map out the energy levels in these delicate quantum systems.

One important technique is photoluminescence excitation (PLE) spectroscopy, which measures the tiny optical signals produced by single-photon emitters when they are excited by continuous-wave (cw) laser light.

### Wide frequency range

“Researchers typically want to measure the response from the sample over a wide frequency range,” explains Jaroslav Sperling, a chemical physicist at Hübner Photonics. “You need a light source that generates light of a very well defined frequency, and that can easily be tuned across a wide range of frequencies.”

## Company

### HÜBNER Photonics

Through continuous technology development, customer orientation and an ISO certified quality management system, HÜBNER Photonics has become a preferred supplier of lasers to major instrument manufacturers and leading research labs for cutting-edge applications in the areas of fluorescence microscopy, flow cytometry, Raman spectroscopy, metrology, holography, nanophotonics and quantum research. HÜBNER Photonics has manufacturing sites in Kassel, Germany and Stockholm, Sweden with direct sales and service offices in USA and UK.

[www.hubner-photonics.com](http://www.hubner-photonics.com)



The C-WAVE from Hübner Photonics combines wide tunability with a narrow linewidth needed and a higher spectral purity than equivalent tunable lasers.

Over the last few years, however, improved design techniques have yielded more sophisticated and more efficient nonlinear materials, such as periodically-poled lithium niobate, that can be phase-matched to the pump laser. A new generation of continuous-wave lasers also provides higher pump powers across the frequency spectrum, making it possible for the first time to produce widely tunable cw OPOs in a turnkey system that delivers narrow-linewidth output at powers levels of a few hundred milliwatts [1].

### Big advantage

One of the big advantages of OPO technology, compared to a conventional cw tunable laser, is that it provides more convenient control of the output wavelength. “There is no need to change the lasing medium or the laser cavity optics to achieve the desired wavelength, since in principle you can generate almost any colour in the visible range with a carefully chosen and suitably operated non-linear crystal,” explains Sperling.

OPO technology also makes it easier to measure the spectra from different samples with the same experimental set-up. “There is no need to change or realign any of the optical components, because the position and direction of the beam always stays the same,” explains Niklas Waasem, a laser physicist at Hübner Photonics who specializes in quantum applications.

“In contrast, other tunable lasers typically need re-alignment of the beam path or even different optics inside the laser after larger changes in the wavelength,” he continues. “With an OPO, researchers can scan through the full frequency range from a computer, which makes it much quicker to characterize different colour centers in the same sample, as well as different samples or material systems.”

The latest commercial systems – such as Hübner Photonics’ C-WAVE – combine this wide tunability with the narrow linewidth needed to resolve specific spectral resonances. OPOs also offer a higher spectral purity than equivalent tunable lasers, which ensures that the weak optical signals generated by atomic-scale defects are not obscured by amplified spontaneous light emission that lasers can generate at unwanted wavelengths. “The spectral purity of an OPO is a huge advantage for PLE experiments,” says Waasem.

Quantum researchers have been quick to recognize the benefits of continuous-wave OPOs for assessing the potential of different single-photon emitters. For example, scientists from the USA, Germany, China, Singapore, and Japan have exploited the C-WAVE platform to measure the photoluminescence spectra from silicon, germanium

and tin vacancy centers in diamond, none of which suffer from the same sensitivity to electric field as nitrogen defects.

In room temperature experiments carried out on an ensemble of germanium vacancy centers (see graph), a broad sweep through the visible frequency range revealed a strong peak in the PLE spectrum at around 602 nm [2]. More detailed experiments at 5 K explored the excitation dynamics of this specific single-photon emitter, showing that resonant fluorescence is only seen when the vacancy center is also excited by a separate laser that can effectively act as an on/off switch [3].

The C-WAVE system has also been used by researchers in Germany to study individual defect centers in hBN, which are difficult to characterize because they have a wide range of transition energies across the entire visible spec-

## What happens next?

### Intensity modulation takes the next step


Once an ensemble of colour centers has been characterized using photoluminescence spectroscopy using a widely tunable laser light source, researchers often turn to other optical techniques to probe the dynamics of the most promising candidates at a specific frequency. These experiments typically require a light source operating at a single well-defined wavelength, and which also enables the intensity of the light to be changed extremely rapidly.

Until recently the most common approach would be to combine a laser with an acousto-optical modulator. Now, however, the same performance can be achieved with modulated diode lasers, which makes the setup more compact, more robust and easier to align. “Modulated diode lasers from Hübner Photonics are popular for ini-

tializing and reading out spin-states because they offer fast modulation capabilities, a high on/off extinction ratio, a high level of spectral purity, as well as an excellent Gaussian beam profile,” comments Niklas Waasem, laser physicist at Hübner Photonics.



Cobolt 06 MLD laser

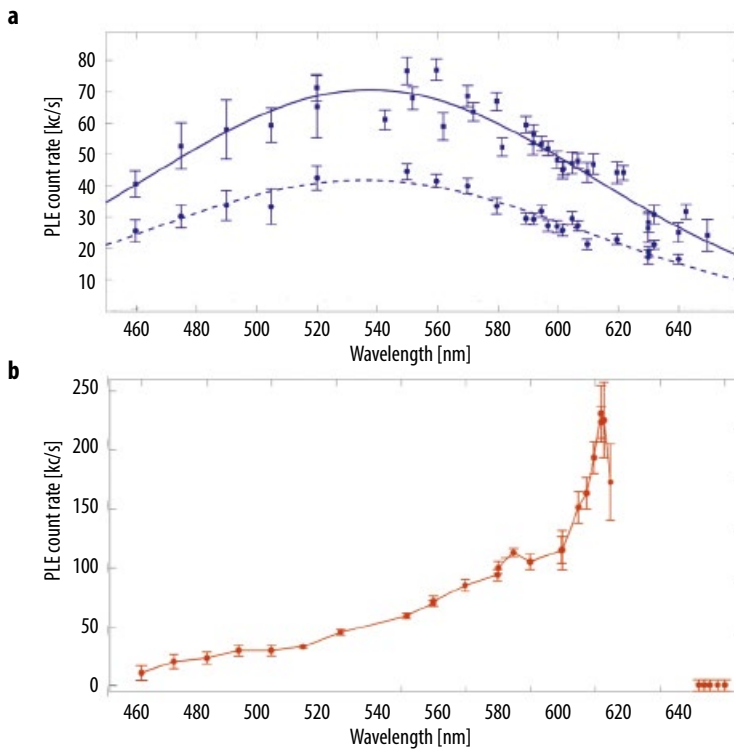

Photoluminescence excitation spectra of an ensemble of Si-V centers recorded at two excitation powers (top). Both spectra are recorded at room temperature and for excitation wavelength tuning from 460 nm to 650 nm (a). Photoluminescence excitation spectrum of an ensemble of Ge-V. The spectrum is recorded at room temperature for excitation wavelength tuning from 450 nm to 640 nm [2] (b).

trum [4]. Measuring the photoluminescence intensity at more than fifty different excitation frequencies revealed the effect of phonon coupling on light emission, which offers a route to identifying the most efficient single-photon emitters in this 2D material.

Such early successes suggest that continuous-wave OPOs will become a popular choice for characterizing a range of different quantum systems. "The OPO concept can be used across different tuning ranges, which will allow for continuous adaption to new experimental requirements," says Sperling. "Overall, we expect continuous-wave OPO technology to mature into a recognized choice among laser light sources that advance the rapidly evolving field of quantum research."

\*

This article was written by Physics World on behalf of Hübner Photonics

- [1] J. Sperling, and K. Hens: Made easy: CW laser light widely tunable across the visible, *Optik & Photonik* 13 (2018) 3, 22-24; DOI: 10.1002/opph.201800022
- [2] S. Häußler et al.: Photoluminescence excitation spectroscopy of SiV- and GeV- color center in diamond, *New J. Phys.* 19 (2017) 6, 063036

- [3] D. Chen et al.: Optical gating of resonance fluorescence from a single germanium vacancy color center in diamond, *Phys. Rev. Lett.* 123 (2019) 3, 033602
- [4] D. Wigger et al.: Phonon-assisted emission and absorption of individual color centers in hexagonal boron nitride, *2D Mat.* 6 (2019) 3, 035006

Contact: Dr Elizabeth Illy  
Director of Marketing

**Hübner Photonics**

phone: +46 8 545 912 30

fax: +46 8 545 912 31

e-mail: [marketing@coboltilasers.com](mailto:marketing@coboltilasers.com)

Web: [www.hubner-photonics.com](http://www.hubner-photonics.com)