


Cobolt 05-01 Series

High Power | Single Frequency | CW Diode pumped lasers

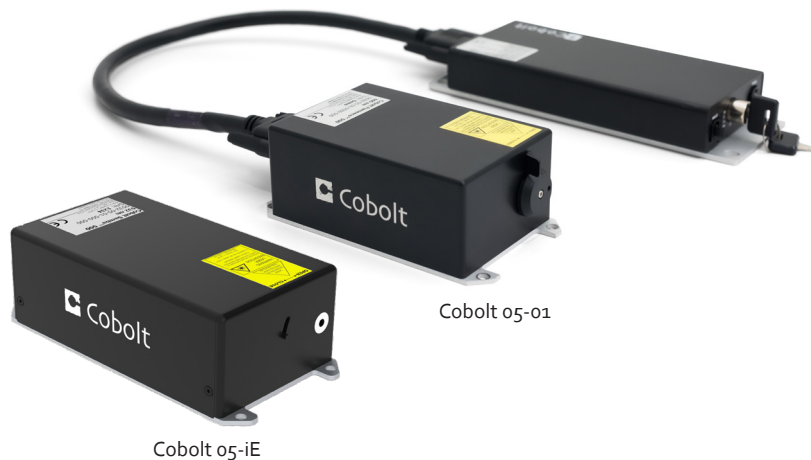

Applications

Raman Spectroscopy
Interferometry
Holography
Particle Analysis
Semiconductor Inspection
Flow Cytometry
Super-resolution Microscopy
Optical trapping
Brillouin Scattering




- CW output power up to 3 W in a perfect beam
- Extremely high level of spectral purity and stability
- Stable single frequency operation over wide temperature range
- Ultra-low intensity noise, down to $< 0.1\%$
- 320 nm, 349 nm, 355 nm, 457 nm, 473 nm, 491 nm, 515 nm, 532 nm, 561 nm, 640 nm, 660 nm, 785 nm, 830 nm and 1064 nm
- Fully Integrated electronics option available
- Up to 24 months warranty, unlimited hours

The Cobolt 05-01 Series lasers are continuous-wave diode pumped laser (DPL) devices operating at a fixed wavelength between 320 nm and 1064 nm. The lasers are built using proprietary HTCure™ manufacturing technology for ultra-robustness in a compact hermetically sealed package.

The Cobolt 05-iE is a fully integrated laser device, including all control electronics. The Cobolt 05-iE eliminates the need for an external controller, bringing the trusted laser performance of Cobolt 05-01 Series into a compact, self-contained device.

The lasers emit a very high-quality laser beam with stable characteristics over a wide range of operating conditions. Single frequency operation provides a narrow spectral bandwidth and long coherence length. The lasers are designed and manufactured to ensure a high level of reliability.

The Cobolt 05-01 Series lasers are intended for stand-alone use in laboratory environments or for integration as OEM components in instruments for applications including fluorescence microscopy, flow cytometry, DNA sequencing, semiconductor inspection, Raman spectroscopy, interferometry, holography and particle analysis.

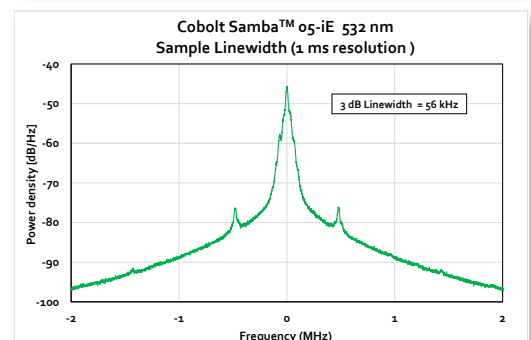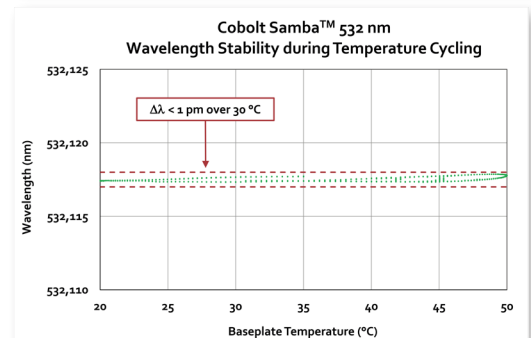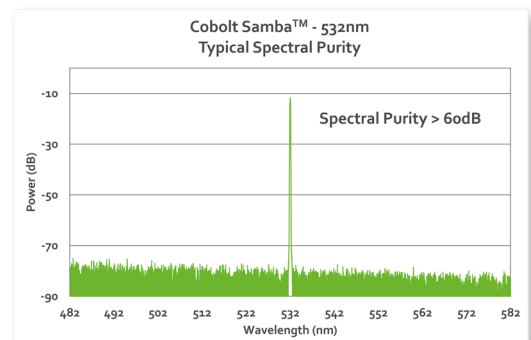

HÜBNER Photonics



Cobolt 05-01 Series

Performance Specifications

	Zydeco™ *	Kizomba™ *	Zouk™ *	Twist™	Blues™	Calypso™	Fandango™
Wavelength in air (nm)	319.8 ± 0.3	348.9 ± 0.3	354.8 ± 0.3	457.0 ± 0.3	473.0 ± 0.3	491.5 ± 0.3	514.8 ± 0.3
Power (mW) without isolator [with intergrated isolator] (mW)	20 [n/a]	50 [n/a]	10 20 [n/a]	100 200 300 [n/a]	100 200 300 [n/a]	200 [n/a]	300 [n/a]
Power stability (±2°C and 8hrs)	< 2%						
Noise, 20 Hz - 20 MHz (pk-pk)	< 7 %		< 2 %			< 5 %	< 2 %
Noise, 20 Hz - 20 MHz (rms)	< 0.7 %		< 0.2 %			< 0.5 %	< 0.2 %
Beam diameter at aperture (µm)	700 ± 50						
Beam symmetry at aperture	> 0.90:1						
Beam divergence (full angle, mrad)	< 0.8			< 1.1			< 1.2
Spatial mode (TEM ₀₀)	M² < 1.1						
Spectral linewidth (FWHM)	< 500 kHz						
Wavelength stability (±2°C and 8hrs)	< 1 pm						
Polarization ratio (linear, vertical)	> 100:1						
Warranty (unlimited hours)	12 months					24 months	12 months

	Samba™	Jive™	Bolero™	Flamenco™	Disco™ *		Rumba™
Wavelength in air (nm)	532.1 ± 0.3	561.2 ± 0.3	639.6 ± 0.3	659.6 ± 0.3	785.0 ± 0.3	830.0 ± 0.3	1064.2 ± 0.6
Power (mW) without isolator [with intergrated isolator] (mW)*****	500 1000 1500** 1800*** [n/a]	300 500 750 1000*** [n/a]	300 500 [n/a]	300 500 [n/a]	500 [400]	500 [n/a]	500 **** 1000 **** 2000 ***** 3000 [n/a]
Power control (% of nominal power)	[n/a]				1 - 100 %		[n/a]
Power stability (±2°C and 8hrs)	< 2%						
Noise, 20 Hz - 20 MHz (pk-pk)	< 1%		< 7 %	< 1%	< 3 %		< 1%
Noise, 20 Hz - 20 MHz (rms)	< 0.1%		< 1 %	< 0.1%	< 0.3 %		< 0.1%
Beam diameter at aperture (µm)	700 ± 50 1000 ± 50**	700 ± 50			1000 ± 100		1000 ± 50
Beam symmetry at aperture	> 0.95:1						
Beam divergence (full angle, mrad)	< 1.2		< 1.4	< 1.5	< 1.3		< 1.6
Spatial mode (TEM ₀₀)	M² < 1.1						M² < 1.2
Spectral linewidth (FWHM)	< 500 kHz				< 100 kHz		< 500 kHz
Wavelength stability (±2°C and 8hrs)	< 1 pm						
Polarization ratio (linear, vertical)	> 100:1						
Warranty (unlimited hours)	24 months***	24 months***	12 months	24 months	12 months		24 months

* Cobolt Zydeco™ 320 nm, Kizomba™ 349 nm, Zouk™ 355 nm and Disco™ lasers available only as 05-01.
** Cobolt Samba™ 1500 mW is available with either a 700 µm or 1000 µm beam dimeter. Model configuration -X01.
*** Cobolt Samba™ 1800 mW and Jive™ 1000 mW warranty is limited to 12 months.
**** Fiber pigtailed option available, contact a sales representative for more information.
***** For more information on the Disco with integrated optical isolator contact a sales representative.

Communication Interface

Communication	USB or RS-232
Standard Baudrate	115200

Model Number

WWWW-05-XX-PPPP-CCCC

Wavelength ↑ ↑ Power ↑ Configuration:

Version: _____

01 Free beam, no optical isolator
03 Fiber pigtailed, no optical isolator****
11 Free beam, integrated optical isolator*****

Configuration:

500 = Gen 5b Controller, RS-232, CE / CDRH
600 = Gen 5b Controller, RS-232, OEM
700 = Gen 5b Controller, USB, CE / CDRH
800 = Gen 5b Controller, USB, OEM
1100 = Integrated electronics, CE / CDRH
1200 = Integrated electronics, OEM
XX01 = 1 mm beam**
XXXX = OEM customization



Fiber pigtailed 05-iE



WARNING VISIBLE AND INVISIBLE LASER RADIATION!



Avoid exposure to beam.
Class 3B Laser Product
Classified per IEC 60825-1:2014

Wvl (nm)	Max.Pwr (mW)
320	200
349	200
355	60
457	499
473	499
491	499
515	499
561	499
660	499



Avoid eye or skin exposure to direct or scattered radiation.
Class 4 Laser Product
Classified per IEC 60825-1:2014

Wvl (nm)	Max.Pwr (mW)
532	3500
561	3000
640	1500
660	1000
785	2000
830	2000
1064	4000

Cobolt 05-01 Series

Operational Environment


The optical performance specifications are not effected by the choice of electronics configuration. However when choosing between the Cobolt 05-01 and 05-iE the operation environment, power supply requirements and thermal management must be considered.

	05-01	Zydeco, Kizomba Bolero	05-iE
Standard power supply*	15 VDC, 6 A		12 VDC, 6.7 A
Voltage acceptance range	11 - 28 V	15 - 28 V	11.2 - 13.2 V
System power consumption	< 65 W, typical 30W		
Maximum laser head baseplate temperature	50 °C	45 °C	45 °C
Ambient temperature, operation	10 - 40 °C	10 - 35 °C	10 - 35 °C
Laser head heatsink thermal impedance (at max ambient temperature)	< 0.2 K/W	< 0.18 K/W	< 0.15 K/W
Beam pointing stability (over operation temperature range)	< 10 µrad/°C, typical 5 µrad/°C		
Ambient temperature, storage	-10 -> +60 °C		
Humidity	0 - 60 % RH non-condensing		
Ambient air pressure	950 - 1050 mbar		



* Cobolt lasers are tested and delivered with the standard power supply.

Electrical Interfaces

Cobolt 05-01 - Laser head

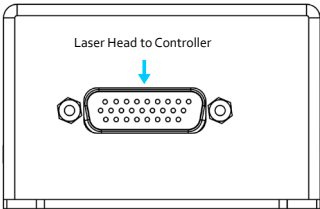


Cobolt 05-01 - Controller

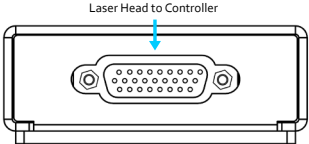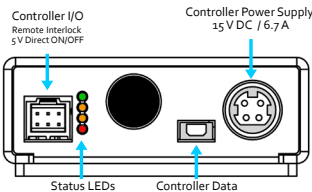

Molex 6 pin - Controller I/O

Pin	Function
1	Remote interlock
2	0 V – Ground
3	Direct Input
4	--
5	LED 1 (LASER ON)
6	LED 2 (ERROR)

Cobolt 05-iE - Laser head

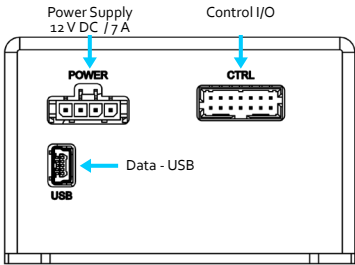


Molex 14 pin- Control I/O

Pin	Function
1	Remote interlock
2	0 V – Ground
3	0 V – Ground
4	RS-232 TX
5	RS-232 RX
6	LED 1A (LASER ON)
7	LED 1B (LASER ON)
8	LED 2 (ERROR)
9	--
10	--
11	Key Switch
12	Direct Input
13	0 V – Ground
14	--

Molex 4 pin - Power Supply

Pin	Function
1	0 V – Ground
2	0 V – Ground
3	+ 12 V - DC
4	+ 12 V - DC

Cobolt 05-iE - Key control box

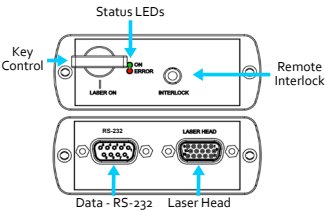

Sub-D 15 pin- Control I/O

Pin	Function
1	LED 1A (LASER ON)
2	LED 2 (ERROR)
3	--
4	0 V – Ground
5	Key Switch
6	--
7	RS-232 TX
8	RS-232 RX
9	--
10	0 V – Ground
11	Remote interlock
12	--
13	--
14	--
15	0 V – Ground


Sub-D pin- RS-232

Pin	Function
1	--
2	RS-232 TX
3	RS-232 RX
4	--
5	0 V – Ground
6	--
7	--
8	--
9	--


Cobolt 05-01 Series

Mechanical Specifications


Cobolt 05-01 Laser head

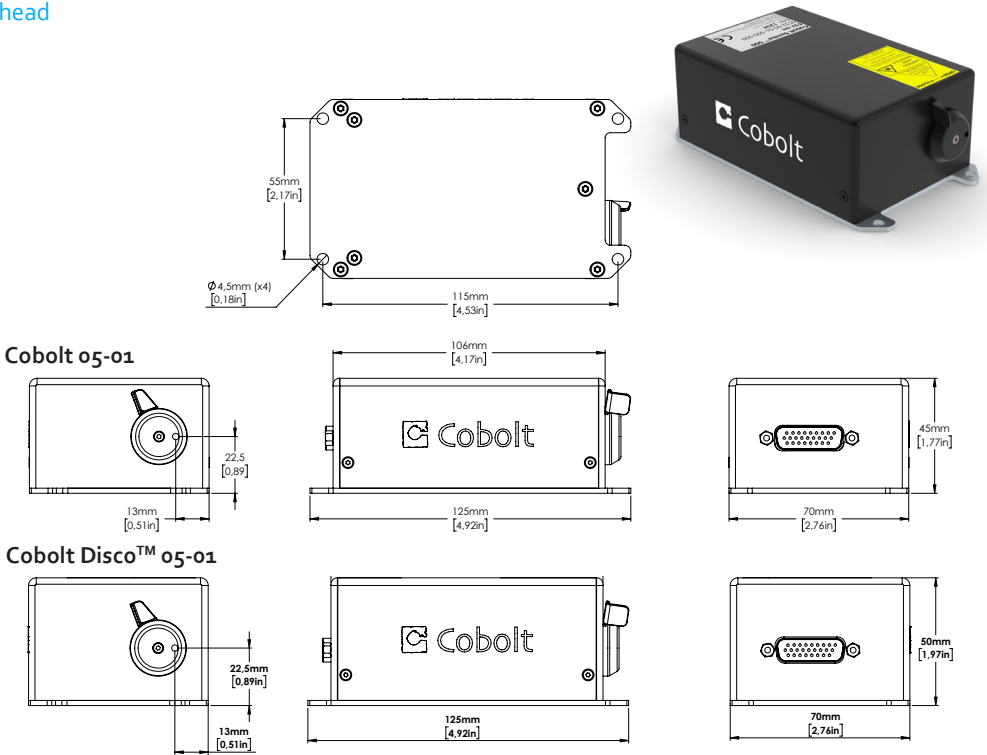

Cobolt 05-11 Disco Laser head with integrated optical isolator

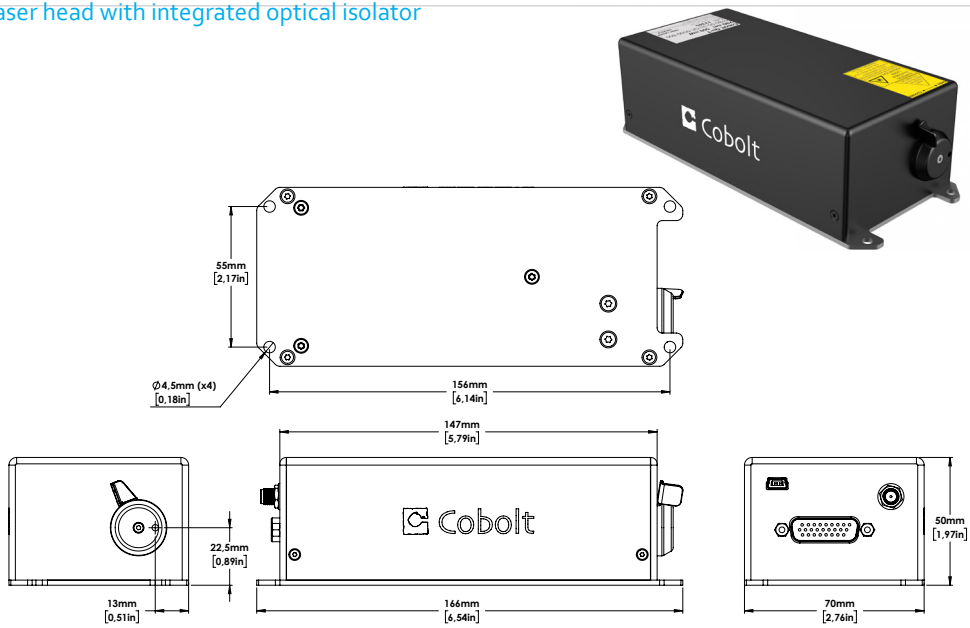


Cobolt 05-01 - Controller

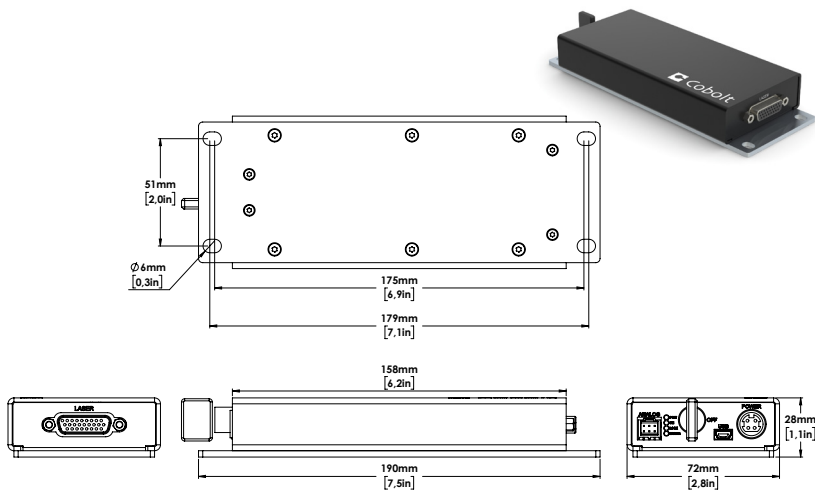


Cobolt 05-01 Series

Mechanical Specifications

Cobolt 05-iE Laser head

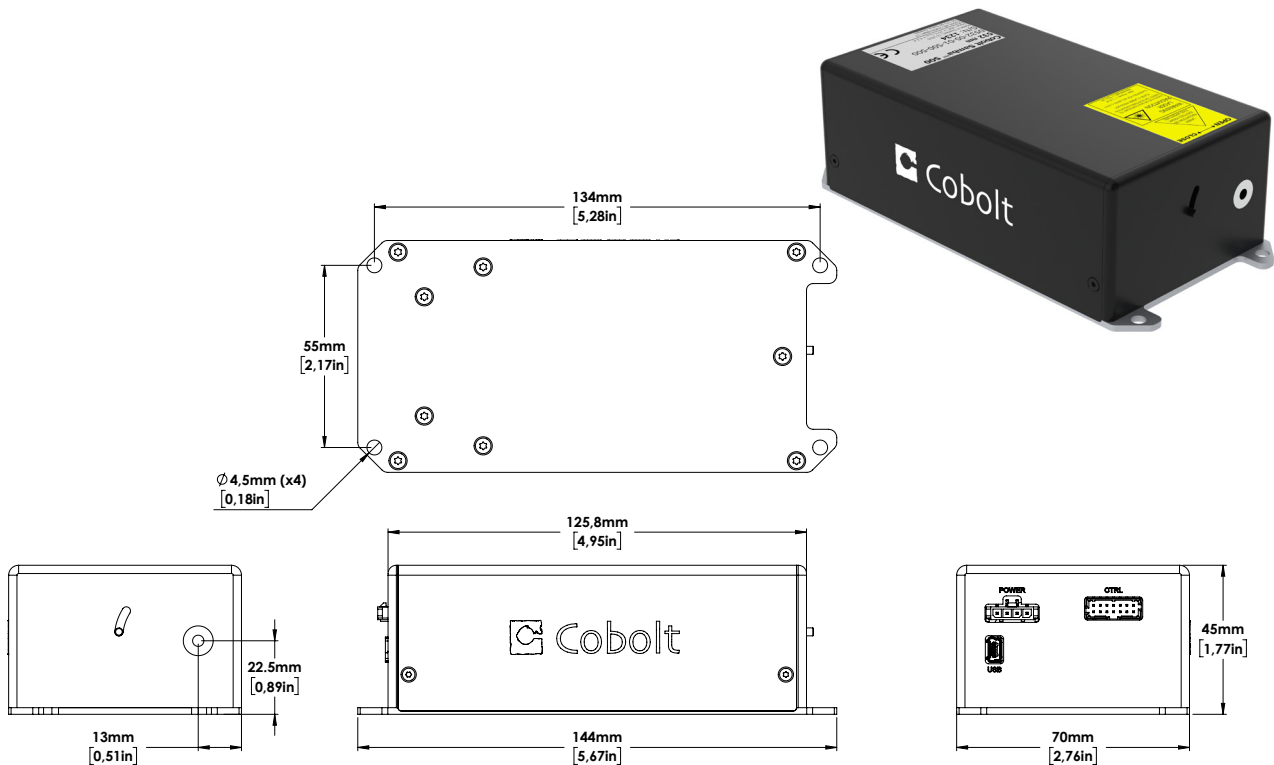

Cobolt 05-iE - Key control box

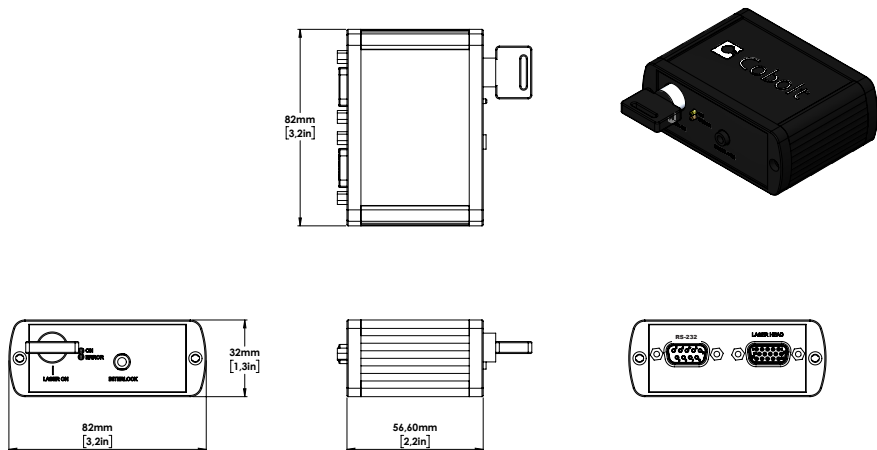

Cobolt 05-01 Series

Options and Accessories


- C-FLEX Laser combiner
- Laser head heatsink with fans for 05-01 lasers : HS-04
- Laser head heatsink with fans for 05-iE lasers : HS-05
- TEC Plate for active baseplate temperature control
- Heatsink with fiber coupling for 05-01 lasers : FIC-04




C-FLEX Laser combiner



Heatsink with fans

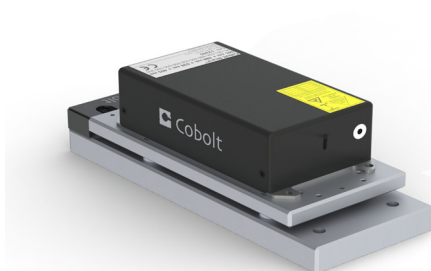

TEC-Plate for active baseplate temperature control

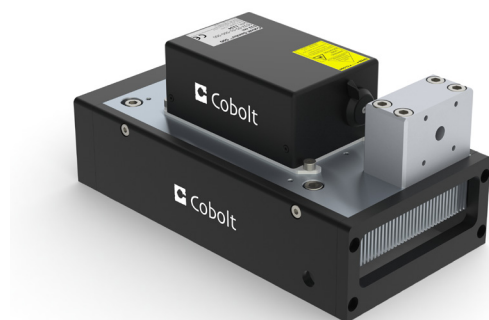

Heatsink with fans for fiber coupling FIC-04

Our Locations

Cobolt AB, a part of HÜBNER Photonics
(Sales in Norway, Sweden, Finland and Denmark)
Solna, Sweden
Phone: +46 8 545 912 30
Fax: +46 8 545 912 31
E-mail: info.se@hubner-photonics.com

HÜBNER Photonics GmbH
(Sales in Germany, Switzerland and Austria)
Kassel, Germany
Phone: +49 561 994 060-0
Fax: +49 561 994 060-13
E-mail: info.de@hubner-photonics.com

HÜBNER Photonics Inc.
(Sales in USA, Canada and Mexico)
San Jose, California, USA
Phone: +1 (408) 708 4351
Fax: +1 (408) 490 2774
E-mail: info.usa@hubner-photonics.com

HA Photonics Pty Ltd
(Sales in UK and Ireland)
London
United Kingdom
Phone: +44 7359 440 871
E-mail: info.uk@hubner-photonics.com

VALO Innovations, a part of HÜBNER Photonics
(VALO Sales and Service)
Hannover, Germany
Phone: +49 511 260 390 70
E-mail: info.valo@hubner-photonics.com

[Find local sales representatives at hubner-photonics.com](http://hubner-photonics.com)